

DNA Analyst Training Laboratory Training Manual

Subject 4: DNA Amplification



PRESIDENT'S
DNA
INITIATIVE



Table of Contents

Purpose.....	3
Objectives	3
Preparation for Exercises	3
Literature.....	4
Exercise 1: Thermal Cyclers.....	5
Purpose.....	5
Tasks	5
Resources	5
Exercise 2: Amplification of Dilution Series.....	6
Purpose.....	6
Tasks	6
Resources	6
Exercise 3: Amplification of Two Donor Mixed Samples	7
Purpose.....	7
Tasks	7
Resources	7
Exercise 4: Amplification of Previously Extracted Samples.....	8
Purpose.....	8
Tasks	8
Resources	8
Subject Review	9
Optional Exercise 5: Effects of Annealing Temperature on Amplification	10
Purpose.....	10
Tasks	10
Resources	10
Exercise Review Questions.....	10
Optional Exercise 6: Effects of Increasing Extension Time on Amplification	11
Purpose.....	11
Tasks	11
Resources	11
Exercise Review Questions.....	11
Optional Exercise 7: Amplification of Degraded Samples.....	12
Purpose.....	12
Tasks	12
Resources	12
Exercise Review Questions.....	12
Optional Exercise 8: Amplification of Inhibited Samples.....	13
Purpose.....	13
Tasks	13
Resources	13
Exercise Review Questions.....	13

Purpose

To instruct the trainee on the amplification process, including the proper techniques for amplification set-up and thermal cycler operation and maintenance.

The following subjects are interdependent:

- Subject 4: DNA Amplification
- Subject 5: Amplified DNA Product Separation
- Subject 6: STR Data Analysis and Interpretation

Upon the sequential completion of these subject areas, the trainee will be able to interpret the data produced in these laboratory exercises.

The trainee will also be introduced to the laboratory's procedures for quality control, to include procedures for limiting the risk of contamination.

Objectives

Upon successful completion of these exercises, the trainee will be able to:

- Describe the theories and procedures of the amplification process.
- Describe the controls used in the amplification process.
- Describe the reagents and their functions in the amplification process.
- Demonstrate the operation of the thermal cycler, including maintenance procedures.
- Explain the conditions under which mineral oil is used in the amplification process.
- Describe the quality system procedures employed by the laboratory to avoid, detect, and document contamination.
- Perform the laboratory's amplification process, per laboratory SOPs.

Preparation for Exercises

Trainer Responsibilities

1. Provide documented safety practices specific to chemicals used in the amplification process, including the pertinent MSDSs.
2. Assign required samples to be amplified as outlined in the Individual Training Plan.
Note: Ensure that the trainee saves samples from each exercise for subsequent exercises.
3. Demonstrate each amplification procedure, including modifications.
4. Demonstrate the proper use of a thermal cycler.
5. Demonstrate and discuss the proper calibration and maintenance of a thermal cycler.
6. Observe the trainee perform the proper calibration and maintenance of a thermal cycler.
7. Observe the trainee performing each step in the amplification procedure.
8. Determine the assessment criteria.
9. Review, verify, and document exercise completion.

Trainee Responsibilities

1. Review documented safety practices specific to chemicals used in the amplification process, to include pertinent MSDSs.
2. Observe each step of the amplification procedure(s), including modifications.
3. Perform each amplification exercise.
4. Perform the calibration and maintenance of a thermal cycler.
5. Document and submit exercise completion, as required by the trainer.

Literature

[Return to Laboratory Training Manual User Guide](#)

Exercise 1: Thermal Cyclers

Purpose

To observe and perform the proper maintenance, calibration, and operation of each type of thermal cycler per the laboratory's SOP's. This demonstration should include the establishment of cycling parameters, operation, and proper documentation of the maintenance performed on the thermal cycler(s). (Proper documentation may include logbooks and/or worksheets.)

Tasks

- Identify and explain the functions of the various components of the instrument
- Establish and verify the cycling parameters
- Calibrate the thermal cycler, as per laboratory SOPs
- Demonstrate proper operation of the thermal cycler(s)

Resources

Sample Protocols: [4.01](#), [4.02](#)

User Manuals: [Applied Biosystems](#)

Exercise 2: Amplification of Dilution Series

Purpose

To familiarize the trainee with the amplification process and assess the trainee's skills. Following the laboratory's SOPs (including controls and blanks), perform the laboratory's amplification method on a dilution series from a known sample.

Tasks

Using the table below, make dilutions of a single source known sample and amplify:

Dilution Factor	Concentration in ng/μl	Amplification volume in μl	Total Target of DNA in ng
Neat sample	0.5	10	5
1:1 of 0.5 ng/μl	0.25	10	2.5
1:1 of 0.25 ng/μl	0.125	10	1.25
1:1 of 0.125 ng/μl	0.0625	10	0.625
1:1 of 0.0625 ng/μl	0.03125	10	0.3125
1:1 of 0.03125 ng/μl	0.015625	10	0.15625
1:1 of 0.015625 ng/μl	0.0078125	10	0.078125
1:1 of 0.0078125 ng/μl	0.0039062	10	0.039062

Note: Values noted above show the dilutions; the use of significant figures was not taken into account.

Resources

Sample Protocols: [4.01](#), [4.02](#)

User Manuals: [Applied Biosystems](#)

Exercise 3: Amplification of Two Donor Mixed Samples

Purpose

To amplify and collect data from two donor mixed samples for use in Subject 6: STR Data Analysis and Interpretation. This data will be used to determine the percent contribution necessary to detect a minor contributor in two donor mixture samples.

Tasks

Using the table below, make dilutions from two single source known samples and amplify:

Sample Number	Donor #1 (ng)	Donor #2 (ng)
1	.1	.9
2	.3	.7
3	.5	.5
4	.7	.3
5	.9	.1

Resources

Sample Protocols: [4.01](#), [4.02](#)

User Manuals: [Applied Biosystems](#)

Exercise 4: Amplification of Previously Extracted Samples

Purpose

To perform the laboratory's amplification procedure(s) on all samples (including the non-probative, mock case samples) extracted in the extraction section of the Laboratory Training Manual, following the laboratory's SOPs (including controls and blanks).

Tasks

Amplify the following:

- All samples extracted in the extraction section of the Laboratory Training Manual

Resources

Sample Protocols: [4.01](#), [4.02](#)

User Manuals: [Applied Biosystems](#)

Subject Review

After completion of the laboratory manual exercises and having previously completed the corresponding theory modules, the trainee should be able to answer the following questions:

- How is contamination limited during amplification set-up?
- What is the purpose of controls, to include positive, negative, and reagent blanks?
- What are the components of each amplification kit? What is the purpose or function of each component?
- Why do most commercial kits utilize Taq Gold versus Taq?
- What are the steps in the PCR process? What occurs during each step?
- What is the expected effect of decreasing the annealing temperature? Of increasing the annealing temperature?
- What is the expected effect on decreasing the extension time? Of increasing the extension time?
- What is multiplexing?
- How does multiplexing make DNA analysis more efficient?
- What factors must be taken into consideration when designing a multiplex PCR reaction?
- What is the specific chemistry of the multiplexing kit used?
- What is the effect of a mutation in a primer binding site?
- What is the laboratory's current target concentration for amplification of single source samples?
- What is the expected result at the higher DNA template concentration? At the lower DNA template concentration?
- What is preferential amplification?
- What is differential amplification?
- What are the conditions under which differential and/or preferential amplification occurs?
- What inhibitors of amplification may co-extract with DNA.
- How do inhibitors affect amplification?
- What effect, if any, do the extraction reagents have on amplification, to include Chelex® 100 and phenol/chloroform/isoamyl alcohol?
- What procedures can be used to overcome inhibition?
- What is the effect of increasing the template DNA concentration to be amplified (e.g. 5 ng or more)?
- Is DNA amplification performed differently when a mixture is known to be present? Explain.
- What is the expected result when a minor component of a mixture is less than 10% of the total concentration?
- Are there any benefits to using commercial extraction kits?
- What amplification reagents are considered critical reagents per the laboratory's SOPS?
- How is it determined that a thermal cycler is functioning properly during a run?
- What maintenance is regularly performed on a thermal cycler?

Optional Exercise 5: Effects of Annealing Temperature on Amplification

Purpose

To collect data from samples amplified using increased and decreased annealing temperatures for use in Subject 6: STR Data Analysis and Interpretation. This data will be used to determine the effects of increasing and decreasing the annealing temperature (by 10°C).

Tasks

- Amplify several samples using a decreased annealing temperature (10°C below the default).
- Amplify several samples using an increased annealing temperature (10°C below the default).

Resources

Sample Protocols: [4.01](#), [4.02](#)

User Manuals: [Applied Biosystems:](#)

Exercise Review Questions

Upon completing this exercise, the trainee should be able to answer these questions:

- What is the expected effect of decreasing the annealing temperature? Of increasing the annealing temperature?

Optional Exercise 6: Effects of Increasing Extension Time on Amplification

Purpose

To collect data from samples amplified using an increased and decreased final extension time of for use in Subject 6: STR Data Analysis and Interpretation. This data will be used to determine the effects of varying the final extension time.

Tasks

Amplify several samples using increased and decreased final extension times.

Resources

Sample Protocols: [4.01](#), [4.02](#)

User Manuals: [Applied Biosystems](#)

Exercise Review Questions

Upon completing this exercise, the trainee should be able to answer these questions:

- What is the expected effect on decreasing the extension time? Of increasing the extension time?

Optional Exercise 7: Amplification of Degraded Samples

Purpose

To collect data from samples treated with DNase for use in Subject 6: STR Data Analysis and Interpretation. This data will be used to determine the effects on amplification of samples treated with DNase.

Tasks

Using the table below, treat a single source known sample as follows:

Sample Number	Sample Quantity	DNase I Concentration	Time
1	1-2 ng	0.1 unit/ μ l	30 seconds
2	1-2 ng	0.1 unit/ μ l	1 minute
3	1-2 ng	0.1 unit/ μ l	1.5 minutes
4	1-2 ng	0.1 unit/ μ l	2 minutes
5	1-2 ng	0.1 unit/ μ l	5 minutes
6	1-2 ng	0.1 unit/ μ l	20 minutes

Resources

Sample Protocols: [4.01](#), [4.02](#)

User Manuals: [Applied Biosystems](#)

Exercise Review Questions

Upon completing this exercise, the trainee should be able to answer these questions:

- What is the expected effect of the DNase treatment on the extracted DNA? Will this be observed in the amplified product? How?
- What mechanisms are involved with DNA degradation?
- What typing results can be expected from degraded DNA?

Laboratory Training Manual

Subject 4: DNA Amplification

Optional Exercise 8: Amplification of Inhibited Samples

Purpose

To collect data from inhibited samples for use in Subject 6: STR Data Analysis and Interpretation. This data will be used to determine the effects on amplification of inhibited samples.

Tasks

Amplify several samples where the following known inhibitors are added to known or positive control samples:

- 10µl of 10% Chelex® 100
- 10µl of phenol/ chloroform/ isoamyl alcohol

Resources

Sample Protocols: [4.01](#), [4.02](#)

User Manuals: [Applied Biosystems:](#)

Exercise Review Questions

Upon completing this exercise, the trainee should be able to answer these questions:

- What inhibitors of amplification may co-extract with DNA.
- How do inhibitors affect amplification?
- What effect, if any, do the extraction reagents have on amplification, to include Chelex® 100 and phenol/chloroform/isoamyl alcohol?
- What procedures can be used to overcome inhibition?

[Return to Laboratory Training Manual User Guide](#)